



Accessing User Data from BMIF Servers Using FTP:

1. Download a FTP **Client** software (e.g. FileZilla- <https://filezilla-project.org/>)
2. Login to the BMIF website
3. Copy the login details found under 'Access Data' on the Dashboard

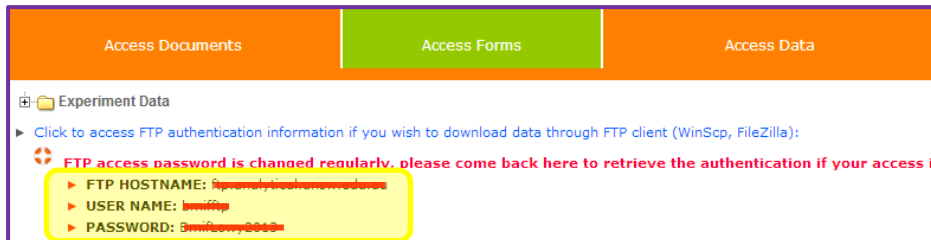


4. Enter the login details into FileZilla and click 'Quick Connect'
5. Navigate to the microscope folder:
 'bmifimages → (Microscope Folder) → (Your data folder)'
6. Drag and Drop folders from the server to your computer

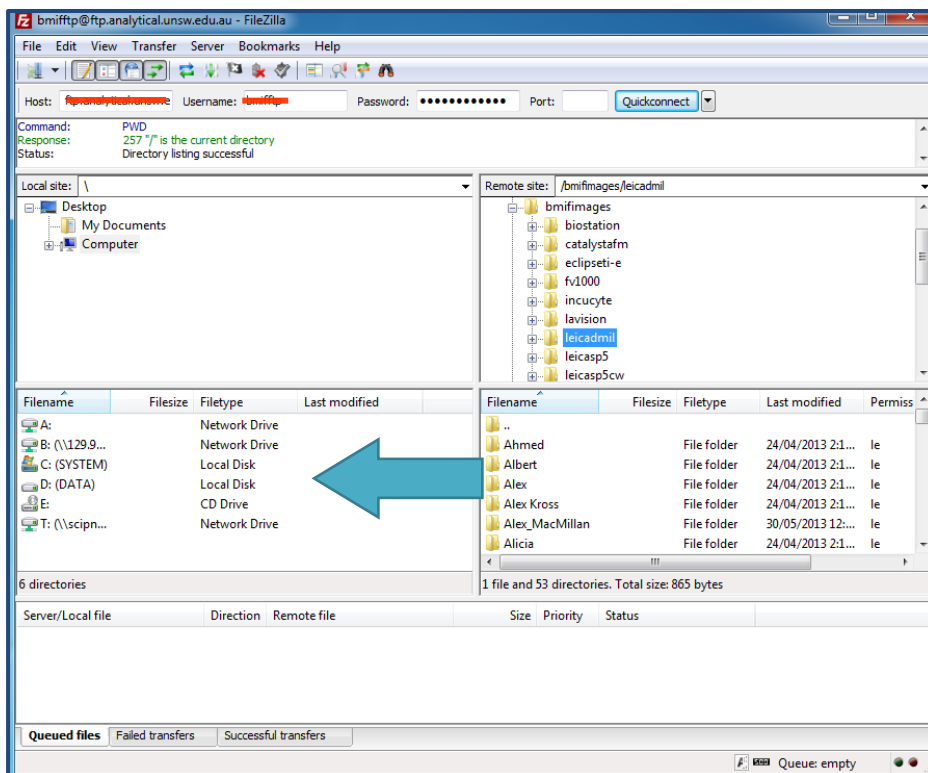

Accessing User Data from BMIF Servers:

1. Login to the BMIF website
2. Select "Access Data" on the Dashboard

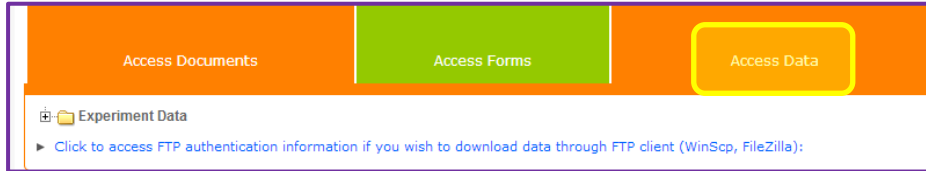

3. Navigate to the microscope folder:
'Experiment Data' → 'FACILITY' → (Microscope Folder)

Experiment Data

- ARCHIVE
- PUBLIC
- FACILITY
 - AFM_Catalyst (Lowy)
 - AFM_MM8 (F10)
 - Biostation
 - Eclipse Ti-E
 - Lavision Intravital
 - Leica DMIL
 - Leica SP5 2P STED
 - Leica SP5 CW STED
 - Nuance
 - Olympus FV1000
 - TILL photonics TIRF
 - Zeiss Axio Examiner
 - Zeiss LSM 780
 - Zeiss PALM
 - Zeiss SD/TIRF

Index of /bmifimages/leicadmil

Name	Size	Date Modified
[parent directory]		
Ahmed/		4/24/13 4:16:00 AM
Albert/		4/24/13 4:16:00 AM
Alex/		4/24/13 4:16:00 AM
Alex_Kross/		4/24/13 4:16:00 AM
Alex_MacMillan/		5/30/13 2:17:00 AM
Alicia/		4/24/13 4:16:00 AM
Anders/		4/24/13 4:16:00 AM
Anjali/		6/20/13 5:29:00 AM
Antonio/		1/6/11 12:00:00 AM
Astrid/		1/4/11 12:00:00 AM
Bakul/		4/24/13 4:16:00 AM
Bin Guan/		4/24/13 4:16:00 AM
Callie/		4/24/13 4:16:00 AM
Caroline/		4/24/13 4:16:00 AM
DMIL Objective Calibrations.txt	865 B	6/11/13 4:42:00 AM
Dacre/		4/24/13 4:16:00 AM
Debie and Peter/		4/18/13 11:42:00 PM
Ellein/		4/24/13 4:16:00 AM
EricQ/		4/24/13 4:16:00 AM
Eve/		4/24/13 4:16:00 AM
Francis/		4/24/13 4:16:00 AM
Gayathri/		4/24/13 4:16:00 AM
Hazim/		4/24/13 4:16:00 AM
Hien/		8/9/13 6:26:00 AM
Hongyu Lu/		4/24/13 4:16:00 AM
Iiveta/		4/24/13 4:16:00 AM
JB/		4/24/13 4:16:00 AM
Jair/		4/24/13 4:16:00 AM
Jarred/		4/24/13 4:16:00 AM
Jenny/		4/24/13 4:16:00 AM
Jessie/Chen/		4/24/13 4:16:00 AM