

# BRIL Staging

MRI

check local regular user uid and gid to make sure the mount matches.

```
$ id nmrsu  
uid=1000(nmrsu) gid=1000(nmrsu) groups=1000(nmrsu),18(dialout),14(uucp),1001(nmruser)
```

**local data located at /data is being synced to:**

```
mkdir /home/nmrsu/ids_bril_staging
```

**service account cred is saved here.**

```
vim /root/.idssvc
```

**For auto mount when startup**

```
vim /etc/fstab
```

to add:

```
//ids.mvac.unsw.edu.au/ids/staging/BRIL/BRIL_MRI /home/nmrsu/ids_bril_staging cifs  
vers=3.0,credentials=/root/.idssvc,uid=1000,gid=1000,forceuid,forcegid,nofail 0 0
```

```
vim /etc/lsyncd.conf
```

to add:

```
# cat /etc/lsyncd.conf  
----  
-- User configuration file for lsyncd.  
--  
-- Simple example for default rsync, but executing moves through on the target.  
--  
-- For more examples, see /usr/share/doc/lsyncd*/examples/  
--  
-- sync{default.rsyncssh, source="/var/www/html", host="localhost", targetdir="/tmp/htmlcopy"}
```

```
settings {
    logfile      = "/var/log/lsyncd.log",
    statusFile   = "/var/log/lsyncd-status.log",
    statusInterval = 20,
}
```

```
sync {
    default.direct,
    source = "/data",
    target = "/home/nmrsu/ids_bril_staging",
    delay  = 10,
    delete = false,
}
```

```
df -h
```

```
//ids.mwac.unsw.edu.au/ids/staging/BRIL/BRIL_MRI 499T 60T 440T 12% /home/nmrsu/ids_bril_staging
```

```
drwxr-xr-x. 2 nmrsu nmrsu 0 Nov 16 16:02 ids_bril_staging
```